



Antibody Stabilizer TRIS

Product description

pH-value: pH 7.2 ± 0.2

Preservative: contains 0.1% ProClin® 300

Storage: shelf life at 2 – 8 °C: 24 months – **Does not tolerate freezing!**

For research use only, not for diagnostic use

Available

Package sizes:	50 ml	order number	130 050
	125 ml	order number	130 125
	500 ml	order number	130 500

Instructions for use

Antibody Stabilizer based on TRIS is made for long-term storage of proteins and antibodies at 2 – 8 °C in solution.

Antibody Stabilizer contains additives for conservation of structure of proteins and antibodies in solution. Thus proteins and antibodies are prevented from losing functionality by storage.

Immediately before using the *Antibody Stabilizer* should be mixed thoroughly.
Just dissolve proteins or antibodies with *Antibody Stabilizer* for storage in the refrigerator.

Stability of proteins and antibodies differs significantly from case to case. Stability depends on characteristics and concentration of the used proteins and antibodies. Thus it can't be predicted generally. User has to test it therefore with its own proteins or antibodies.

Antibody Stabilizer is ready-to-use.

For further information please visit www.candor-bioscience.com.

ProClin is a registered trade mark of Rohm and Haas Company.